# **T400-Series Surgical Protocol**

# Dog Renal Artery: Chronic Blood Flow Measurement

### **APPLICATION BASICS**

Site: Renal Artery

Species: Dog
Weight: 15 kg
Duration: Chronic
Vessel Diameter: 2.5 mm

PROBE

Size: 3 mm Connector: 10-pin


Reflector: L with sliding cover

Cable Length: 1 meter

Catalog #: MC-3PSS-LS-WC100-CRS10-GC

FLOWMETER TS420 Perivascular Module

# Flow Ranges Observed

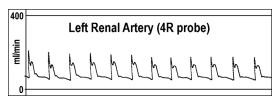

#### **ACKNOWLEDGEMENT**

Dr. Allan Buchholz, Sterling Research Group, 81 Columbia Turnpike,

Rensselaer, NY 12144 Vita Lanoce, Squibb Pharmaceuticals, Rt. 206 and Provinceline

Rd., Princeton NJ. 08543

## **Application**

The measurement of renal blood flow has an important role in research on hemodynamics, electrolyte regulation and pregnancy induced hypertension. This protocol was used to test potential antihypertensive agents. Other studies have focused on diuretics and nephrotoxic agents. While average renal flow may also be obtained from the renal vein, the pulsatile waveform of the renal artery provides additional information and visual confirmation of a functioning chronic implant.

Variations of this protocol include a ventral approach for situations where access to both kidneys is desired or other protocols require a laparotomy. One experienced researcher sutures the Probe with the cable facing the opposite direction and a loose suture suspending the Probe from the body wall. Another reports fixing the position of the kidney by suturing the renal capsule to the body wall.

# **Surgical Approach**

Premedicate with 0.02 mg/kg atropine IM, induce with 18 mg/kg thiamylal and maintain on halothane. With anesthetized dog in right lateral recumbency, palpate the left kidney and visualize the location of the renal vessels. Make a 4 cm vertical skin incision over the kidney. Continue the incision by bluntly dissecting through the most superficial muscle layer, the cutaneous trunci, to expose the external abdominal oblique muscle. The external oblique muscle can easily be identified by the caudoventral direction of its fibers.

Bluntly dissect through it to expose the cranioventral fibers of the internal oblique muscle. Continue the dissection through the internal abdominal oblique and the underlying transversus abdominous muscle. Continue the incision through retroperitoneal fat to expose the kidney and renal vessels. Gently dissect the tissue cranial to the hilus of the kidney and identify the renal vessels. Take care not to damage the ureter exiting caudolaterally. Locate the renal artery and place a silk suture around it. Note that the renal artery often bifurcates. Make sure that there isn't another branch. Place a 3 mm Probe around the artery (or paired arteries) and close the sliding bracket. Securing the Probe

takes special care because the kidney is mobile and suspended by the vascular pedicle. The objective is to fix the position of the

transonic

THE MEASURE OF SETTER RESULTS.


(Continued on next side.)

# Dog Renal Artery: Chronic Blood Flow Measurement Cont.

## Surgical Approach cont.

Probe with respect to the artery without making the Probe support the weight of the kidney. Suture the cable to the body wall close to the point where the renal artery joins the aorta. This position minimizes the strain on the cable when the kidney moves. Place sutures through the bracket suture holes to adjacent connecting tissue. These sutures will limit rotation of the Probe around the artery. Wrapping the Probe with silicone sheet wrap is recommended to stabilize the implant.

Pass the cable through the incision and close each layer of the body wall independently. Create a subcutaneous pouch for the cable and implantable connector. Close the subcutaneous layer and skin routinely. By seven days post surgery, fibrotic tissue effectively seals the tunnel containing the cable. When a flow reading is desired, the dog may be lightly anesthetized to exteriorize the implantable connector.

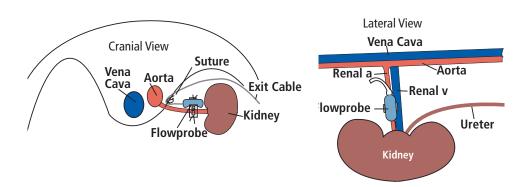

### **REFERENCES**

Mann, W.A., Kinter, L.B., and Woodward, P., "Selective DA-1 Antagonism of SK&F R-87516-J, Fenoldopam, and Dopamine-Stimulated Renal Vasodilation in Dogs", Federation Proceedings, Vol. 45, No. 3, p. 427, 1986.

Mangino, M.J., Jendrisak, M.D., Brunt, E., Anderson, C.B., "Eicosanoid Synthesis Inhibition and Renal Function during Acute Rejection", Transplantation, Vol. 45, No.5, pp. 902-907, 1988.



Transonic Systems Inc. is a global manufacturer of innovative biomedical measurement equipment. Founded in 1983, Transonic sells "gold standard" transit-time ultrasound flowmeters and monitors for surgical, hemodialysis, pediatric critical care, perfusion, interventional radiology and research applications. In addition, Transonic provides pressure and pressure volume systems, laser Doppler flowmeters and telemetry systems.

### **AMERICAS**

Transonic Systems Inc. 34 Dutch Mill Rd Ithaca, NY 14850 U.S.A. Tel: +1 607-257-5300 Fax: +1 607-257-7256

support@transonic.com

### **EUROPE**

Transonic Europe B.V.
Business Park Stein 205
6181 MB Elsloo
The Netherlands
Tel: +31 43-407-7200
Fax: +31 43-407-7201
europe@transonic.com

### ASIA/PACIFIC

Transonic Asia Inc. 6F-3 No 5 Hangsiang Rd Dayuan, Taoyuan County 33747 Taiwan, R.O.C. Tel: +886 3399-5806 Fax: +886 3399-5805 support@transonicasia.com

#### **JAPAN**

Transonic Japan Inc. KS Bldg 201, 735-4 Kita-Akitsu Tokorozawa Saitama 359-0038 Japan Tel: +81 04-2946-8541 Fax: +81 04-2946-8542 info@transonic.jp